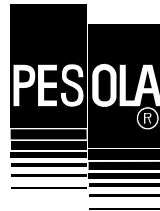
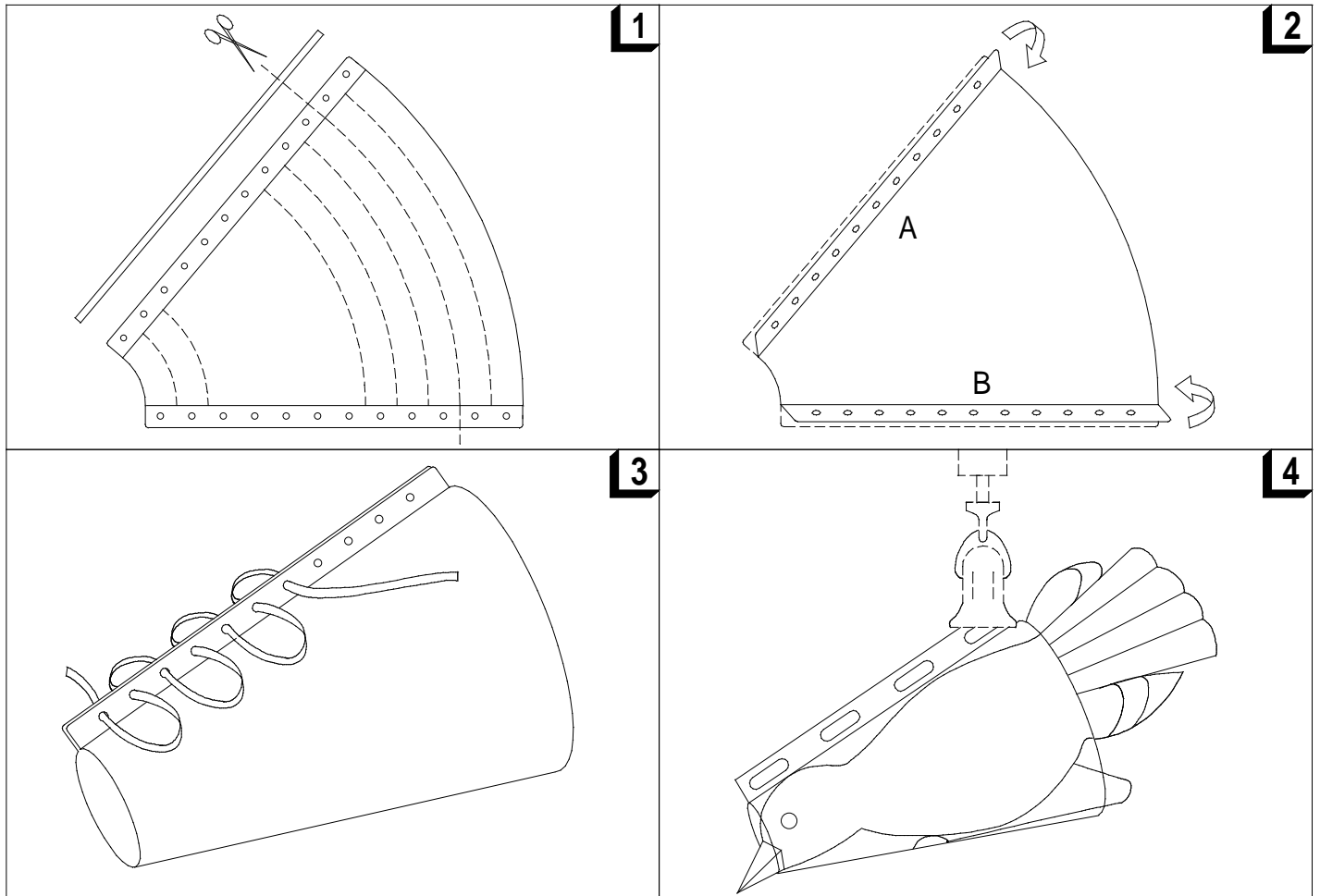
3.655

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## English

### Features

- Suitable for repeated use with 50g - 600 g PESOLA precision scales for field and laboratory application.
- Conical form gently holds wings together and precise measurement is facilitated by immobilization of animal.
- Opening at base of cone allows the animal air and facilitates cleaning.
- Cone can be disassembled after use to save space.
- Also suitable for other small animals.

### Instructions for assembly and use

1. The size of the cone can be reduced by cutting along one of the marked lines to fit very small animals.
2. Fold "A" and "B" at 180° angle and press down firmly.
3. Roll up the foil conically and join opposite holes with nylon tape. Tare the scale in vertical position with cone attached.
4. Place bird head down, legs forward in the cone and with arm supported read the weight. To avoid a falsification of the results protect the scale from wind. Only leave the animal in the cone during the short process of weighing to prevent a build up of heat. Before further weighing remove droppings, check zero position and, if necessary, adjust tare.
5. Do not store cone in direct sunlight.